

## Announcements

---



### Risø TL/OSL Reader software updates available on the web

You may now download the latest software for your TL/OSL reader from the Risø home page [www.osl.risoe.dk](http://www.osl.risoe.dk). Once you reach the website, select the option on the left hand side of the screen for 'The Riso TL/OSL reader' and then the section entitled 'Software'. As well as being able to download the latest release of the different software packages, there is also a section entitled "Revision History" where you can see the most important changes that have been made to the software since previous versions were released.

For instance the Sequence Editor has recently had a few small but very practical changes:

1. When you run a sequence, the grid with the commands does not disappear as it used to, but now it is shown in the lower part of the screen next to the usual communication log window. The cell being executed is marked, and you may freely inspect all the cells, but no changes can be made during the run.
2. In "Sequence Options" you may tick "Sequence file copy" (this is the default setting). When this is done, you automatically save a copy of the sequence file (\*.SEC file) and this is automatically given the same name and path as the data file (\*.BIN file). This file can be opened and used again, but not saved again without changing the name. The idea is that you always store a .SEC file as documentation for how the .BIN file was created. A detailed description of this system is available on the home page under the heading "Tech and Change Notes"

Torben Lapp  
Radiation Research Division, Risø DTU, Denmark

